

### Hydroxycinnamic acids of coffee

**Sample:** coffee, water extract

**Inj. solvent:** 0.5% aq. H<sub>3</sub>PO<sub>4</sub>

**Inj. volume:** 10 µL

**SP:** 50x4.6 Reprosil-Pur Basic C18 (1.9 µm)

**MP:** 10% AcN in 0.5% aq. H<sub>3</sub>PO<sub>4</sub>

**Flow:** 2 mL/min

**Detection:**


— 270 nm

— 330 nm

**Detector cell volume:** 10 µL

**Temperature:** 25 °C

**Pressure:** 210 bar

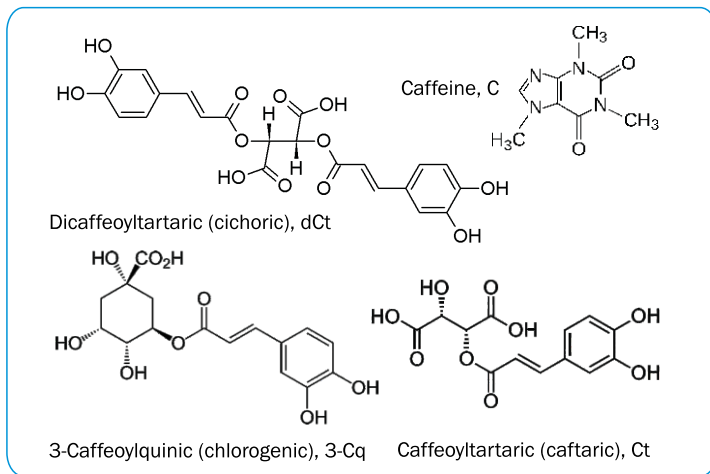

ELSICO HPLC Company



SCAN

Consulting and training for HPLC

[www.hplc.ru/training](http://www.hplc.ru/training)



<https://sites.google.com/site/scansychev/>  
<http://www.chromforum.ru/>