

**Water soluble vitamins**

**Sample:** vitamin containing medicine, sample 2

**Inj. solvent:** mobile phase

**Inj. volume:** 10 uL

**SP:** 150x4.6 Reprosil-Pur C18-AQ (3 um)

**MP:** 5% AcN and 10mM pentasulfonate in 0.1% aq. H3PO4

**Flow:** 2 mL/min

**Detection:**


— 205 nm

— 260 nm

**Detector cell volume:** 10 uL

**Temperature:** 25 °C

**Pressure:** 200 bar

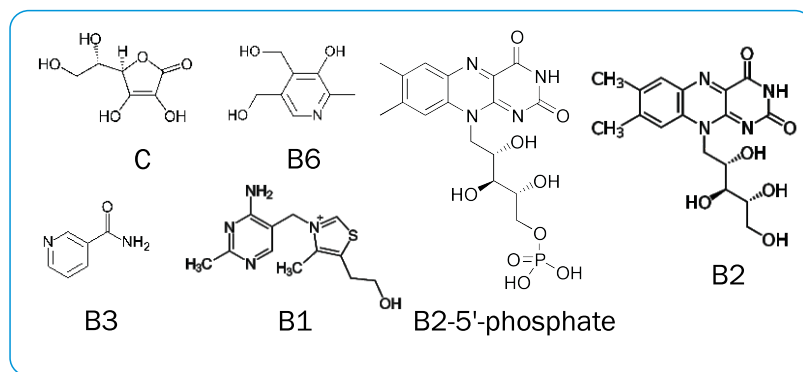

ELSICO HPLC Company



SCAN

Consulting and training for HPLC

[www.hplc.ru/training](http://www.hplc.ru/training)



<https://sites.google.com/site/scansychev/>  
<http://www.chromforum.ru/>